



## AssayMax Human Apolipoprotein C-II ELISA Kit

Catalog Number EA8012-1

### Introduction

Apolipoprotein C-II (ApoC-II) is secreted in plasma and resides in both the very low density lipoproteins (VLDL) and high density lipoproteins (HDL). ApoC-II plays a major role in lipid metabolism as the obligate cofactor for lipoprotein lipase, which catalyzes the hydrolysis of triglyceride-rich lipoproteins (1). The protein has 79 amino acid residues and, in the absence of lipid, self-associates to form amyloid fibrils implicated in the pathogenesis of a number of diseases including Alzheimer's, Parkinson's, and Creutzfeldt-Jakob diseases (2). ApoC-II aggregates are present in human atherosclerotic plaques, and its fibrils initiate macrophage inflammatory responses (3). Deficiency of either apoC-II or lipoprotein lipase results in hypertriglyceridemia (4). Transgenic mice over-expression of human ApoC-II also causes hypertriglyceridemia attributed to the delayed clearance of VLDL triglycerides (5).

### Principal of the Assay

The AssayMax Human Apo C-II ELISA (Enzyme-Linked Immunosorbent Assay) kit is designed for detection of human Apo C-II in plasma, serum, urine, and cell culture samples. This assay employs a quantitative sandwich enzyme immunoassay technique that measures human Apo C-II in less than 4 hours. A polyclonal antibody specific for human Apo C-II has been pre-coated onto a 96-well microplate with removable strips. Apo C-II in standards and samples is sandwiched by the immobilized antibody and biotinylated polyclonal antibody specific for Apo C-II, which is recognized by a streptavidin-peroxidase conjugate. All unbound material is then washed away and a peroxidase enzyme substrate is added. The color development is stopped and the intensity of the color is measured.

### Caution and Warning

- This kit is for research use only.
- The kit should not be used beyond the expiration date.
- The Stop Solution is an acid solution

### Reagents

- **Human Apo C-II Microplate:** A 96-well polystyrene microplate (12 strips of 8 wells) coated with a polyclonal antibody against human Apo C-II.
- **Sealing Tapes:** Each kit contains 3 pre-cut, pressure-sensitive sealing tapes that can be cut to fit the format of the individual assay.
- **Human Apo C-II Standard:** Human Apo C-II in a buffered protein base (1 µg, lyophilized).
- **Biotinylated Apo C-II Antibody (100x):** A 100-fold concentrated biotinylated polyclonal

- antibody against Apo C-II (80 µl).
- **MIX Diluent Concentrate (10x):** A 10-fold concentrated buffered protein base (30 ml).
  - **Wash Buffer Concentrate (20x):** A 20-fold concentrated buffered surfactant (30 ml).
  - **Streptavidin-Peroxidase Conjugate (SP Conjugate):** A 100-fold concentrate (90 µl).
  - **Chromogen Substrate:** A ready-to-use stabilized peroxidase chromogen substrate tetramethylbenzidine (8 ml).
  - **Stop Solution:** A 0.5 N hydrochloric acid to stop the chromogen substrate reaction (12 ml).

## Storage Condition

- Store unopened kit at 2-8<sup>0</sup>C up to expiration date.
- Opened MIX Diluent may be stored for up to 1 month at 2-8<sup>0</sup>C. Store reconstituted reagents at -20<sup>0</sup>C or below.
- Opened unused strip wells may return to the foil pouch with the desiccant pack, reseal along zip-seal. May be stored for up to 1 month in a vacuum desiccator.

## Other Supplies Required

- Microplate reader capable of measuring absorbance at 450 nm
- Pipettes (1-20 µl, 20-200 µl, 200-1000µl and multiple channel)
- Deionized or distilled reagent grade water

## Sample Collection, Preparation and Storage

- **Plasma:** Collect plasma using one-tenth volume of 0.1 M sodium citrate as an anticoagulant. Centrifuge samples at 2000 x g for 10 minutes and assay. Dilute samples 1:500 into MIX Diluent as follows: add 10 µl of sample to 490 µl of MIX Diluent (1:50) to make Solution A; then add 80 µl of Solution A to 720 µl of MIX Diluent (1:10) to make a final working solution (1:500). Store samples at -20<sup>0</sup>C or below for up to 3 months. Avoid repeated freeze-thaw cycles. (EDTA or Heparin can also be used as anticoagulant.)
- **Serum:** Samples should be collected into a serum separator tube. After clot formation, centrifuge samples at 2000 x g for 10 minutes. Dilute samples 1:500 into MIX Diluent as follows: add 10 µl of sample to 490 µl of MIX Diluent (1:50) to make Solution A; then add 80 µl of Solution A to 720 µl of MIX Diluent (1:10) to make a final working solution (1:500). Store samples at -20<sup>0</sup>C or below for up to 3 months. Avoid repeated freeze-thaw cycles.
- **Cell Culture Supernatants:** Centrifuge cell culture media at 2000 x g for 10 minutes to remove debris. Collect supernatants and assay. Store samples at -20<sup>0</sup>C or below. Avoid repeated freeze-thaw cycles.
- **Urine:** Collect urine using sample pot. Centrifuge samples at 800 x g for 10 minutes and assay. Dilute samples 1:4 into MIX Diluent. Store samples at -20<sup>0</sup>C or below for up to 3 months. Avoid repeated freeze-thaw cycles.

## Reagent Preparation

- Freshly dilute all reagents and bring all reagents to room temperature before use. If crystals have formed in the concentrate, mix gently until the crystals have completely dissolved.
- **MIX Diluent Concentrate (10x):** Dilute the MIX Diluent 1:10 with reagent grade water. Store for up to 1 month at 2-8<sup>0</sup>C.

- **Standard Curve:** Reconstitute the 1 µg of Apo C-II Standard with 0.5 ml of MIX Diluent to generate a solution of 2 µg/ml. Allow the standard to sit for 10 minutes with gentle agitation prior to making dilutions. Prepare triplicate standard points by serially diluting the standard solution (2 µg/ml) 1:4 with MIX Diluent to produce 0.5, 0.125, 0.031, 0.078 and 0.0019 µg/ml solutions. MIX Diluent serves as the zero standard (0 µg/ml). Any remaining solution should be frozen at -20°C.

Standard Point	Dilution	[Apo C-II] (µg/ml)
P1	Standard (2 µg/ml)	2.000
P2	1 part P1 + 3 part MIX Diluent	0.500
P3	1 part P2 + 3 part MIX Diluent	0.125
P4	1 part P3 + 3 part MIX Diluent	0.031
P5	1 part P4 + 3 part MIX Diluent	0.0078
P6	1 part P4 + 3 part MIX Diluent	0.0019
P7	MIX Diluent	0.000

- **Biotin Apo C-II Antibody (100x):** Spin down the antibody briefly and dilute the desired amount of the antibody 1:100 with MIX Diluent. Any remaining solution should be frozen at -20°C.
- **Wash Buffer Concentrate (20x):** Dilute the Wash Buffer Concentrate 1:20 with reagent grade water.
- **SP Conjugate (100x):** Spin down the SP Conjugate briefly and dilute the desired amount of the conjugate 1:100 with MIX Diluent. Any remaining solution should be frozen at -20°C.

## Assay Procedure


- Prepare all reagents, working standards and samples as instructed. Bring all reagents to room temperature before use. The assay is performed at room temperature (20-30°C).
- Remove excess microplate strips from the plate frame and return them immediately to the foil pouch with desiccant inside. Reseal the pouch securely to minimize exposure to water vapor and store in a vacuum desiccator.
- Add 50 µl of Apo C-II standard or sample per well. Cover wells with a sealing tape and incubate for two hours. Start the timer after the last sample addition.
- Wash five times with 200 µl of Wash Buffer. Invert the plate and decant the contents, and hit it 4-5 times on absorbent paper towel to completely remove liquid at each step.
- Add 50 µl of Biotinylated Apo C-II Antibody to each well and incubate for one hour.
- Wash five times with 200 µl of Wash Buffer. Invert the plate and decant the contents, and hit it 4-5 times on absorbent paper towel to complete remove liquid at each step.
- Add 50 µl of Streptavidin-Peroxidase Conjugate to each well and incubate for 30 minutes. Turn on the microplate reader and set up the program in advance.
- Wash five times with 200 µl of Wash Buffer.
- Add 50 µl of Chromogen Substrate per well and incubate for about 10 minutes or till the optimal blue color density develops. Gently tap plate to ensure thorough mixing and break the bubbles in the well with pipette tip.
- Add 50 µl of Stop Solution to each well. The color will change from blue to yellow.
- Read the absorbance on a microplate reader at a wavelength of 450 nm **immediately**. Please note that after the reaction is stopped for about 10 minutes, some black particles may be generated at high concentration point, which will reduce the readings.

## Data Analysis

- Calculate the mean value of the triplicate readings for each standard and sample.
- To generate a Standard Curve, plot the graph using the standard concentrations on the x-axis and the corresponding mean 450 nm absorbance on the y-axis. The best-fit line can be determined by regression analysis using log-log or four-parameter logistic curve-fit.
- Determine the unknown sample concentration from the Standard Curve and multiply the value by the dilution factor.

## Standard Curve

- The curve is provided for illustration only. A standard curve should be generated each time the assay is performed.

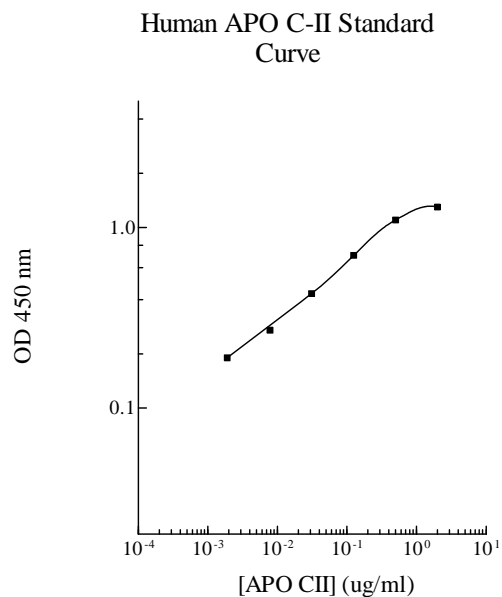

## Performance Characteristics

- The minimum detectable dose of Apo C-II is typically 10 ng/ml.
- Intra-assay and inter-assay coefficients of variation were 4.9% and 7.5% respectively.
- No significant cross reactivity with Apo AI, Apo AII, Apo B, Apo CI, Apo CIII or Apo E.

## Linearity

Sample Dilution	Average Percentage of Expected Value	
	Plasma	Serum
1:250	105%	108%
1:500	100%	97%
1:1000	94%	95%

Sample Dilution	Average Percentage of Expected Value
	Urine
No Dilution	113%
1:2	95%
1:4	92%

## Recovery

Standard Added Value	0.05 – 0.5 ug/ml
Recovery %	82 - 110
Average Recovery %	96

## Cross-Reactivity

Species	% Cross Reactivity
Beagle	< 1
Bovine	None
Monkey	< 10 (suggest 1:10 dilution for plasma)
Mouse	< 2 (suggest 1:4 for plasma)
Rat	< 1
Swine	None

**Plasma reference value: 30 – 80 µg/ml.**

## References

- (1) Jackson CL *et al.* (1984) *Proc. Natl. Acad. Sci. USA* 81:2945-2949
- (2) Hatters, DM *et al.* (2000) *Biochemistry* 39:8276-8283(5)
- (3) Medeiros LA *et al.* (2004) *J. Biol. Chem.* 279:10643-10648
- (4) Fojo SS and Brewer HB (1992) *J. Intern. Med.* 231:669-677
- (5) Shachter NS *et al.* (1994) *J. Clin. Invest.* 93:1683-1690

Version 3.2